


Quick Installation Guide Compass 315M




Complete demo system ships in 2 boxes.

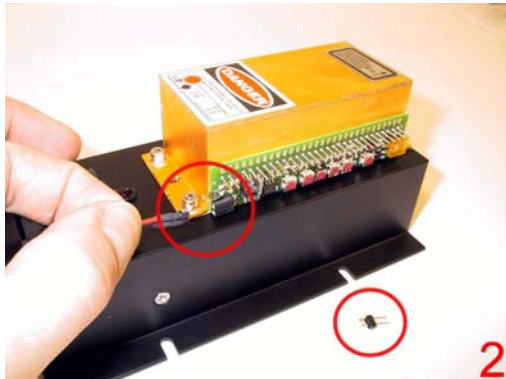

Please use thin layer of thermal compound between laser head and heat sink. Connect fan cable after extra connector (red top, dark bottom) has been removed.



Complete assembled laser head on heat sink.

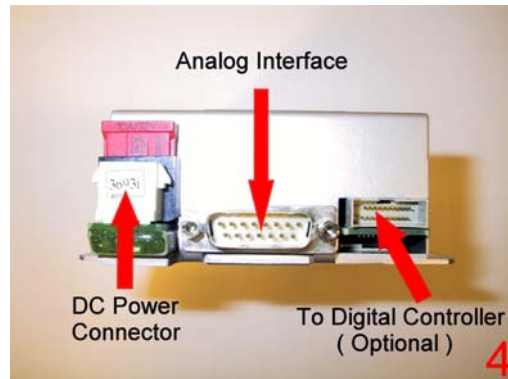

Controller with DC Power Connector, analog interface and connector to digital controller.



Assembled Controller with auto-start adapter, head cable and DC power cable.



Fully assembled system. Please use same ground for laser head and controller. Assemble system before connecting AC or DC Power.



The Compass™ 315M laser is a modular component, sold for use in OEM equipment and is not to be used as a stand-alone laser. The OEM is responsible for compliance with all applicable safety regulations. Please see Compass 315M manual on pages 1-1 to 1-5 for laser safety instructions.

Repackaging Guide Compass 315M




Please attach antistatic foam and plastic cap for protection during transportation.



Arrangement before shipping. (Box 1)



Please put all components back into bags!
(Except head cable and power supply)

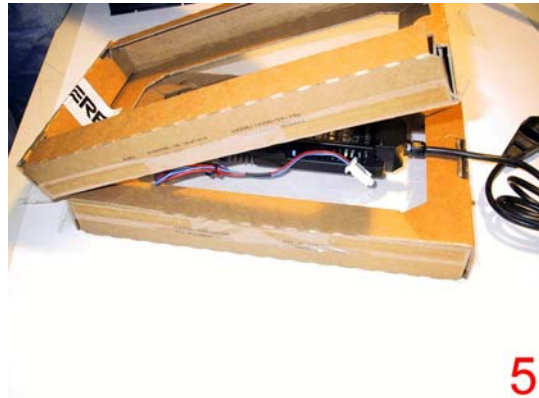

Put power supply and heat sink in Box 2.
Also place devices **between** foils!



Please put head and controller **between** foils for maximum protection in box 1.



Arrangement before shipping. (Box 2)

The Compass™ 315M laser is a modular component, sold for use in OEM equipment and is not to be used as a stand-alone laser. The OEM is responsible for compliance with all applicable safety regulations. Please see Compass 315M manual on pages 1-1 to 1-5 for laser safety instructions.