

Software Installation Guide
ModeMaster PC[™]



Software Installation Guide
ModeMaster PC



27650 SW 95th Ave.
Wilsonville, OR 97070

This document is copyrighted with all rights reserved. Under the copyright laws, this document may not be copied in whole or in part or reproduced in any other media without the express written permission of Coherent, Inc. Permitted copies must carry the same proprietary and copyright notices as were affixed to the original. This exception does not allow copies to be made for others, whether or not sold, but all the material purchased may be sold, given or loaned to another person. Under the law, copying includes translation into another language.

Coherent, the Coherent Logo, and ModeMaster PC are trademarks or registered trademarks of Coherent, Inc. All other trademarks or registered trademarks are the property of their respective owners.

Patents referenced in this document were active as of the printing date of the manual (see last page for date). The patents referenced herein may have expired. You are advised to check to see if the patents are still active: <http://portal.uspto.gov/external/portal/pair>.

Every effort has been made to ensure that the data given in this document is accurate. The information, figures, tables, specifications, and schematics contained herein are subject to change without notice. Coherent makes no warranty or representation, either expressed or implied, with respect to this document. In no event will Coherent be liable for any direct, indirect, special, incidental, or consequential damages resulting from any defects in its documentation.

Technical Support

In the U.S.:

Should you experience difficulties with your product, or need technical information, visit our website: www.Coherent.com. You can obtain additional support by either telephoning our Technical Support Hotline at 1.800.343.4912, or e-mailing our Support Team at support.instruments@Coherent.com. Telephone coverage is available Monday through Friday (except U.S. holidays).

If you call outside our office hours, your call will be taken by our answering system and will be returned when the office reopens.

If there are technical difficulties with your laser that cannot be resolved by support mechanisms outlined above, e-mail, or telephone Coherent Technical Support with a description of the problem and the corrective steps attempted. When communicating with our Technical Support Department via the web or telephone, the model and Laser Head serial number of your laser system will be required by the Support Engineer responding to your request.

Outside the U.S.:

If you are located outside the U.S., visit our website for technical assistance, or telephone our local service representative. Representative phone numbers and addresses can be found on the Coherent website: www.Coherent.com.

Coherent provides web and telephone technical assistance as a service to its customers and assumes no liability thereby for any injury or damage that may occur contemporaneous with such services. These support services do not, under any circumstances, affect the terms of any warranty agreement between Coherent and the buyer. Operating a Coherent product with any of its interlocks defeated is always at the operator's risk.

TABLE OF CONTENTS

| | |
|------------------------------------|----------|
| Preface | v |
| Publication Updates | v |
| Software Installation | 1 |
| Installing the Software | 1 |
| Uninstalling the Software | 2 |

Preface

This guide includes instructions for installing the ModeMaster PC software and USB driver.

For detailed information on ModeMaster PC, refer to the *ModeMaster PC User Manual* (part number 0217-802-00), available in Adobe® Acrobat® PDF format on the installation CD.

Publication Updates

To view information that may have been added or changed since this publication went to print, connect to www.Coherent.com.

ModeMaster PC Software Installation Guide

SOFTWARE INSTALLATION

This section explains how to install or uninstall the ModeMaster PC software.

Installing the Software

Note: The USB driver is automatically installed during this process.

Step 1: Insert the ModeMaster PC CD into your CD-ROM drive.

Step 2: If the installation process does not automatically start, browse to your CD-ROM drive (typically, **D:**) and double-click *Setup.exe*.

Step 3: Follow the on-screen instructions.

*Note: If a window similar to the one shown below appears, click the **Install this driver software anyway** button to continue the installation. (This example is from a Windows 7 installation.)*



Step 4: When the following screen appears, click the **OK** button to complete the software installation.



Uninstalling the Software

To uninstall the software, click the **Start** button in Windows and then click **All Programs > Coherent > ModeMaster PC > Uninstall ModeMaster PC**.

ModeMaster PCTM Software Installation Guide
©Coherent, Inc., 8/2015, printed in the USA
Part No. 1188027 Rev. AB